



Operating Instructions

863 Valley View Road, Eighty Four, PA 15330 USA
Tel: 724-941-9701 • www.skcinc.com

NEED HELP?

SCAN OR CLICK HERE
TO CONTACT US



Air-Met Scientific
www.airmet.com.au

Tedlar® Sample Bags Cat. No. 231-01-TCLP

Cat. No. 231-01-TCLP is a disposable 1-liter, Tedlar film sample bag with a single stainless steel fitting designed for use with a Zero Headspace Extractor (ZHE) for the EPA TCLP method. The use of Tedlar sample bags with stainless steel fittings has been validated by the EPA.

Operation

The bag is designed to be attached directly to the extractor vessel via a special stainless steel coupler.

To open the fitting, remove the knurled screw cap on the fitting. Attach the coupler to the bag by screwing onto the bag fitting, then attach directly to the ZHE unit.

Note: SKC Cat. No. 231-01-TCLP bags are designed for single use applications and are not intended for the storage and transport of hazardous materials.

Tedlar is a registered trade name of DuPont Corporation.

SKC Limited Warranty and Return Policy

SKC products are subject to the SKC Limited Warranty and Return Policy, which provides SKC's sole liability and the buyer's exclusive remedy. To view the complete SKC Limited Warranty and Return Policy, go to <http://www.skcinc.com/warranty>.

www.skcinc.com